


Procedure to confirm upgradable period registered into HASP Key

1. Start HASP Info


It is possible to confirm the upgradable period registered into the HASP key on the screen of Utility software called HASP Info. There are following two ways to start "HASP Info".

(a) Start from the Windows start menu

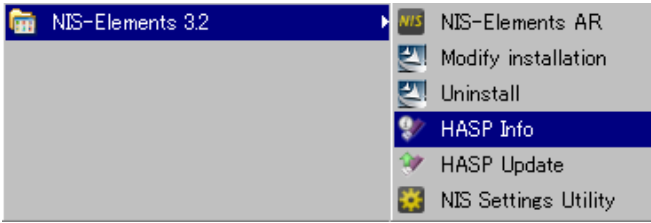

Select "HASP Info" of NIS-Elements from Windows start menu.

(b) Start from NIS-Elements help menu

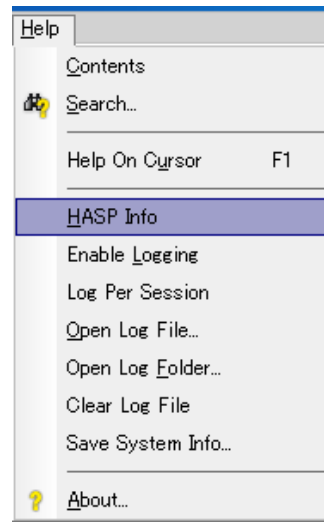

Select "HASP Info" from Help menu of NIS-Elements.

2. Confirmation of the upgradable period

The upgradable period displayed on a HASP Info screen can be confirmed with the following procedures.

HASP ID: 2710771B

Compatible with NIS-Elements 3.2 and higher

HASP 4 memory

		Version
Systems	NIS Elements AR	3.1
	NIS Elements BR	3.1
Modules	ND (6 dimensions)	

HASP SRM memory

		Version
Systems	NIS Elements AR	3.2
	NIS Elements BR	3.2
Modules	ND (6 dimensions)	

Activation Date: August 02, 2010 (1) Activation Date

Upgrades available until September 02, 2011 (2) Upgradable Period

The following information is shown at the bottom of "HASP Info".

- (1) Activation Date
- (2) Upgradable Period

If the expiry date of upgradable period registered into HASP Key is on or after the upgradable term reference date of NIS-Elements^{*1} you can upgrade to the version.

If the expiry date of upgradable period registered into HASP Key is before the upgradable term reference date of NIS-Elements, it is necessary to purchase SUA option for the required period to extend an upgradable period.

*1: You can check the upgradable term reference date of each version of NIS-Elements at the download corner of NIS-Elements. Please refer to next page "3.Confirmation for release date of NIS-Elements" for more detailed information.

3. Confirmation for upgradable term reference date of NIS-Elements

You can check the upgradable term reference date of each version of NIS-Elements at download for NIS-NIS-Elements. In the example of the following figure, the date is displayed in red box is upgradable term reference date of NIS-Elements.

Software	Release date	Download page
Imaging Software NIS-Elements AR Ver4.40.00 for 64bit edition Use Condition 1. Upgrade period - Please check "HASP Info" to confirm if "upgradable term reference date" is April 24, 2015 or later. - Please refer to "The Confirmation Procedure of an upgrade period" for more detailed information. (PDF: [Link]) 2. Supported OS - Microsoft Windows 7 Professional SP1 (English/Japanese). Since V4.00, Windows XP and Windows Vista are not supported. Modification 1. New Function - Supported New Volume View - Supported 3D Object Measurement	01/05/2015	Download (881.5MB)

Summary

The following figures are the summary about the upgradable period registered into the HASP Key, upgradable term reference date, and SUA option.

